



## Microscopes

- ❑ Collect acrylic covers
- ❑ Eyepieces
  - Clean with q-tip or glass cloth and lens cleaner
  - Make sure they're attached firmly
  - Make sure black rubber is in proper position
- ❑ Focus
  - Focus scope
  - Check that focus knobs are securely attached
- ❑ Lenses
  - Rotate to the side (if needed) when replacing specimens
  - Clean lens with q-tip, kimwipe, or glass cloth and lens cleaner if image is foggy
- ❑ Acrylic boxes
  - Keys to open box in info desk top drawer
  - Clean box with soft rag and acrylic cleaner
    - Greasy nose prints
    - Fingerprints around doors

# Microscope Anatomy

Volunteer  
*Academy*

---

